

# Development of a high-throughput CUT&RUN platform for epigenomic mapping of rare primary immune cells


Keith Maier<sup>1</sup>, Ellen N. Weinzapfel<sup>1</sup>, Allison R. Hickman<sup>1</sup>, Matthew R. Marunde<sup>1</sup>, Carolina P. Lin<sup>1</sup>, Danielle N. Maryanski<sup>1</sup>, Liz Albertorio-Saez<sup>1</sup>, Dughan J. Ahimovic<sup>2</sup>, Michael Bale<sup>2</sup>, Juliana J. Lee<sup>3</sup>, Michael-Christopher Keogh<sup>1</sup>, Immunological Genome Project Consortium

<sup>1</sup>EpiCypher Inc., Research Triangle Park, Durham NC 27709, USA. <sup>2</sup>Laboratory of Epigenetics and Immunity, Department of Pathology and Laboratory Medicine, Weill Cornell Medicine, New York, NY, USA. <sup>3</sup>Department of Immunology, Harvard Medical School, Boston, MA, USA.



## Epigenetic regulation is central to cell and gene therapy, but has been challenging to study


- Many genomic strategies for cell & gene therapy focus on transcription; however, RNA-seq reveals the **outcomes** – not driving **mechanisms**
- **Epigenomics is the solution:** Mapping the location of histone post-translational modifications (PTMs) and chromatin-associated proteins, such as transcription factors, provides molecular insights that are central to cell fate and function
- However, existing epigenomic technologies, such as **ChIP-seq**, are limited by high costs, poor sensitivity & reliability, and complicated sample prep
- These challenges have precluded epigenomic analysis for cell & gene therapy

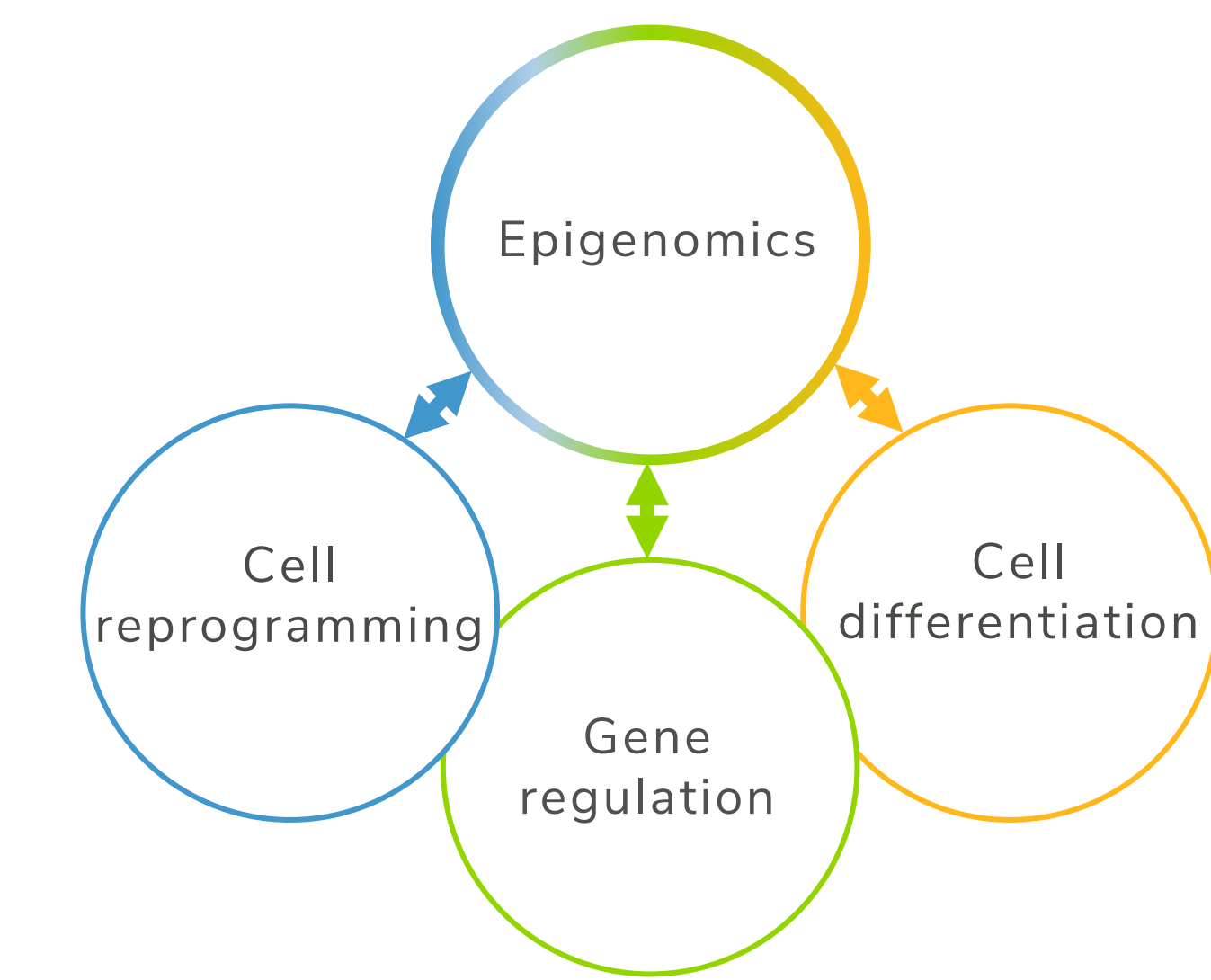

**Figure 1:** Understanding epigenetic regulation is critical to successful cell and gene therapy applications:

- iPSCs
- CAR T-cells
- T cell exhaustion
- dCas9/Cas9 targeting

## CUTANA™ CUT&RUN assays provide important advantages compared to ChIP-seq

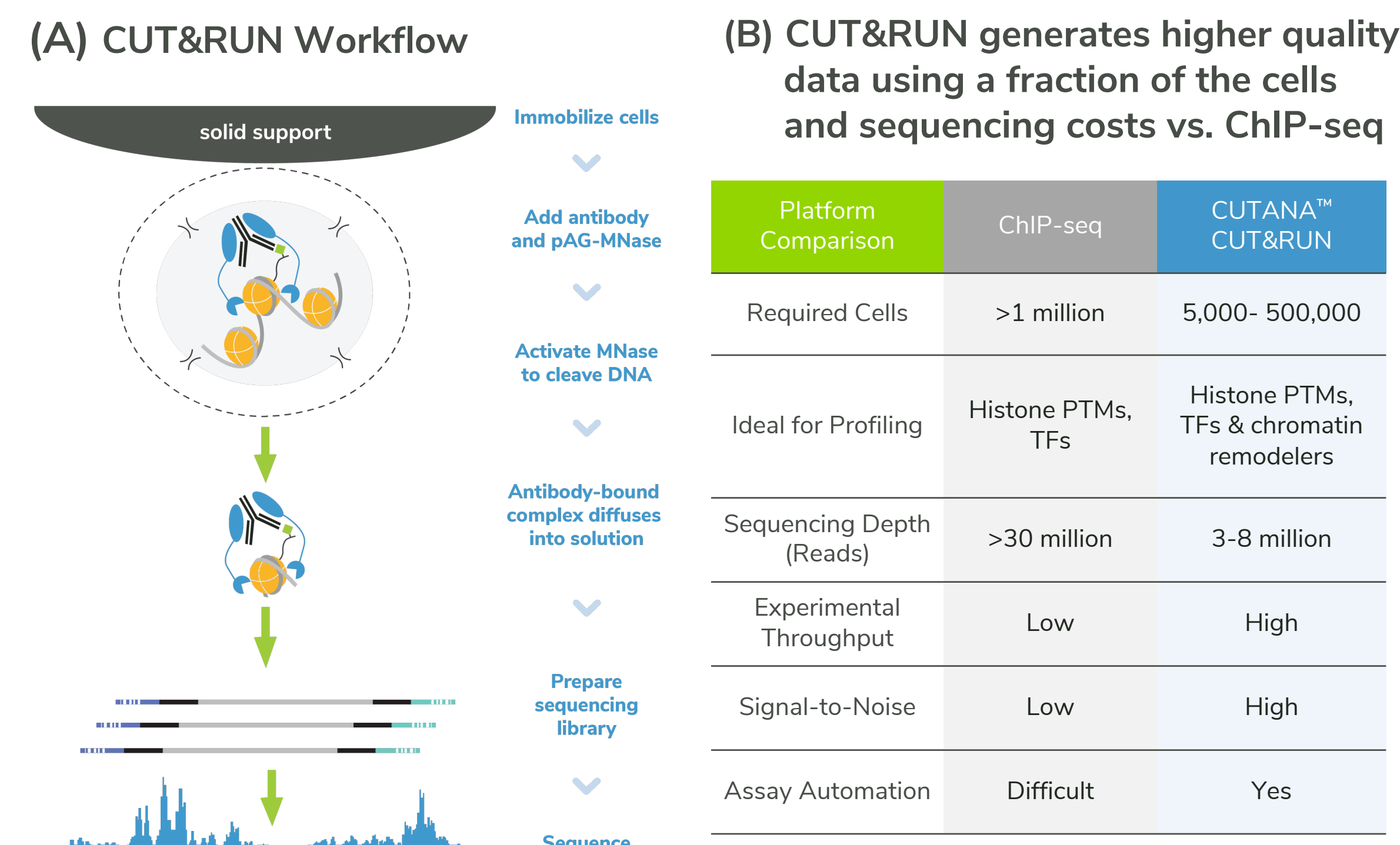

**Figure 2.** Overview of the CUTANA CUT&RUN workflow and advantages compared to ChIP-seq. Because CUT&RUN releases antibody-bound fragments into solution (A), it has improved signal-to-noise even with significantly reduced cell numbers and sequencing depth (B).

## Automated CUTANA™ CUT&RUN Assays: high-throughput, customized epigenomics

**(A) Cut hands-on time in half while increasing throughput 4X**

Freedom EVO100 robotics setup

Platform Comparison	Standard CUT&RUN	Automated CUT&RUN
Length of protocol	4 days	3 days
Total hands-on time	17 hours	8.5 hours
Experimental throughput	24 reactions	96 reactions
Compatible with immune cells?	Yes	Yes
Number primary immune cells	Down to 5,000 (select targets)	Optimized for 10,000

**(B) Processes up to 384 CUT&RUN reactions per week**

### Optimization steps:


- Miniaturized reaction volumes and sequencing library prep for 96-well plates
- Standardized workflow for native & cross-linked cells & nuclei
- Developed spike-in controls to monitor workflows, validate antibodies, and flag failed reactions



Buffer optimization reduces sample loss

## autoCUT&RUN enables ultra-sensitive epigenomic profiling from low cell numbers


**(A) High signal-to-noise down to 5k cells**      **(B) Reliable signal across varying cell numbers**

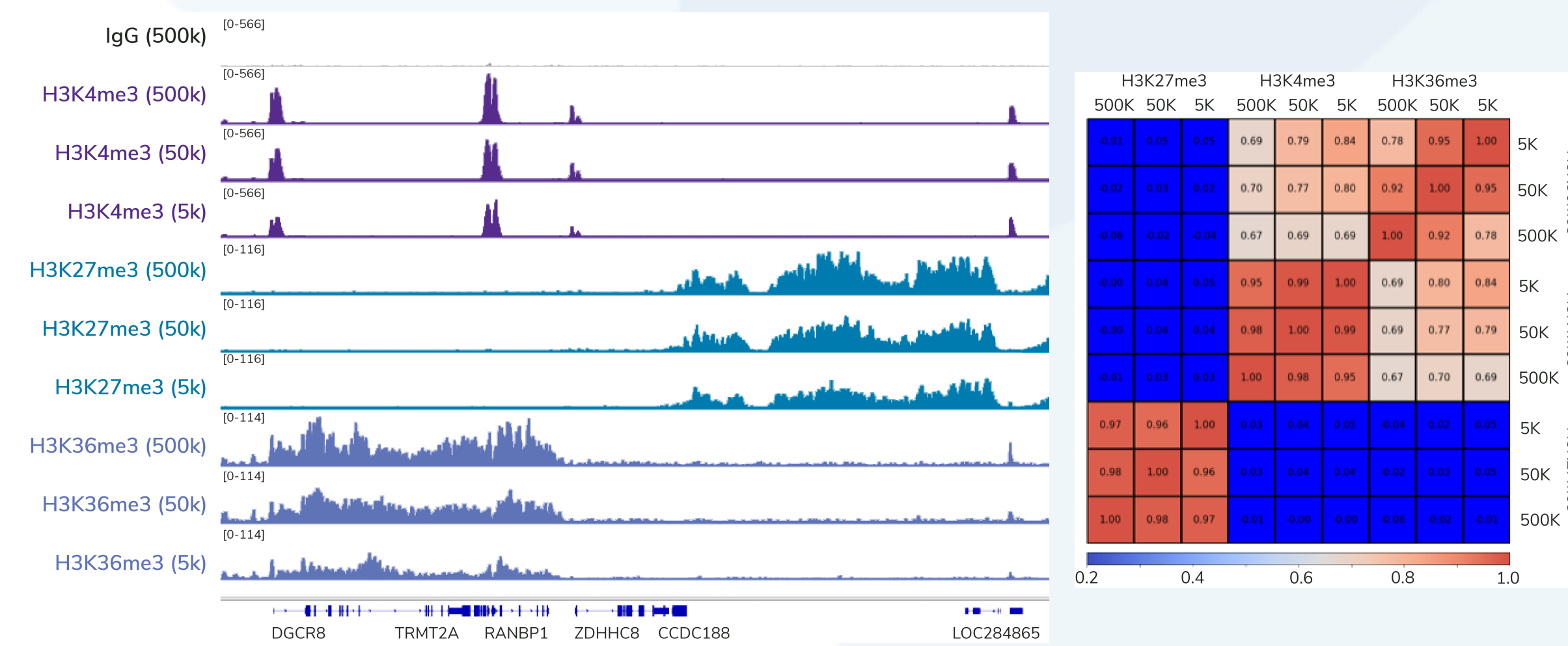

**Figure 4.** Automated CUTANA CUT&RUN (autoCUT&RUN) was used to generate maps for various histone PTMs using decreasing amounts of K562 cells (A). H3K4me3 Pearson correlation matrix (B) shows high concordance across cell numbers for each target.

## The key to autoCUT&RUN: SNAP-CUTANA™ Spike-in Controls and highly specific antibodies


**(A) SNAP-CUTANA Spike-ins for CUT&RUN**

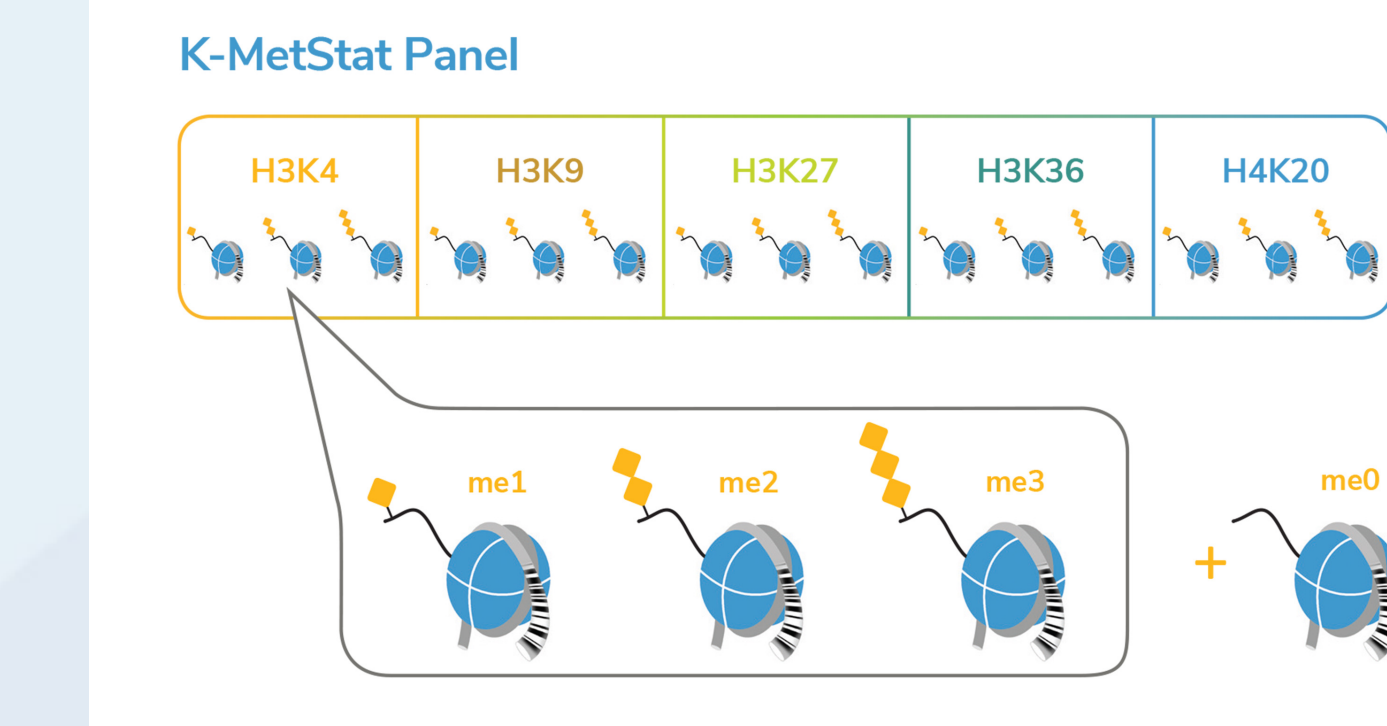

**Figure 5A:** SNAP-CUTANA Spike-ins are panels of DNA-barcoded nucleosomes that are:

- Added to CUT&RUN just prior to antibody addition
- Processed alongside sample as an ideal internal control
- Recovery of on- and off-target spike-ins determines experimental success

**(B) SNAP Spike-ins identify specific histone PTM antibodies for CUT&RUN**

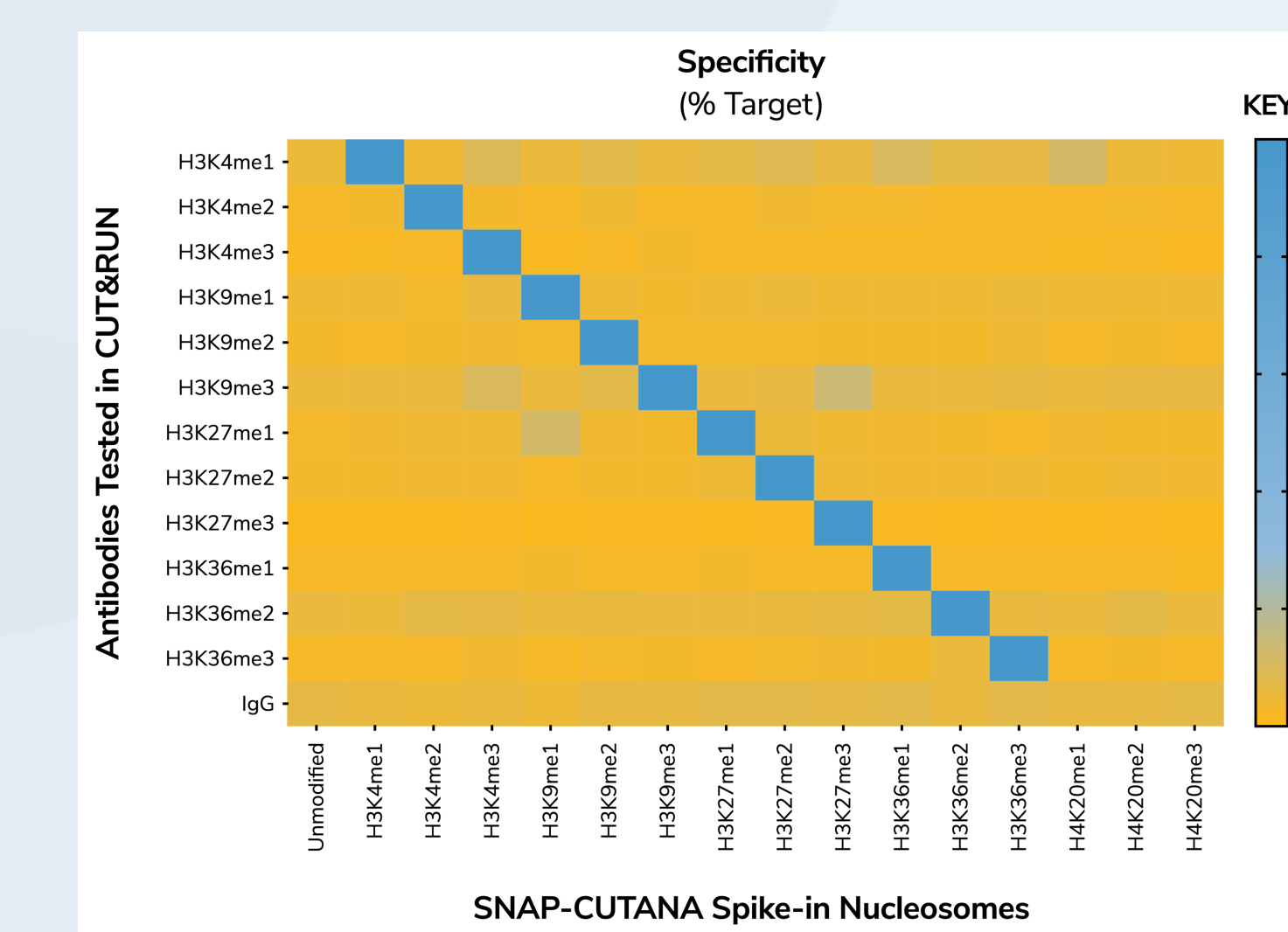


**Figure 5B:** SNAP-CUTANA Spike-ins were used to identify best-in-class CUT&RUN antibodies for widely studied histone lysine methylation marks:

- In the heatmap, each row is an antibody tested in CUT&RUN
- Columns show recovery of individual spike-ins relative to the on-target PTM (yellow indicates low recovery, blue shows cross-reactivity >20%)

## Unique validation strategy enables reliable profiling of transcription factors and chromatin-associated proteins

### Validation of EpiCypher antibody to FOXA1

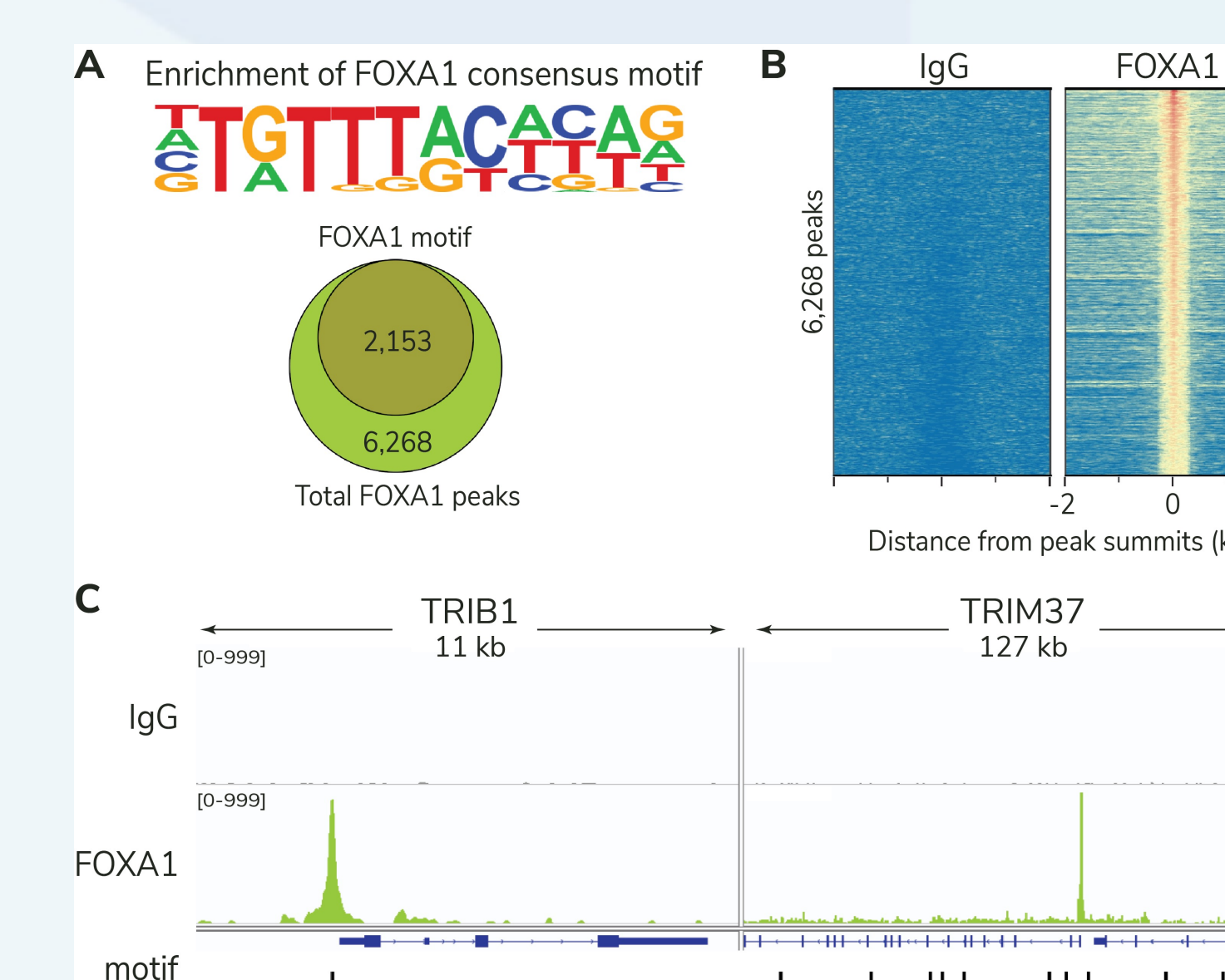


**Figure 6:** Our stringent validation protocol involves multiple genome-wide analysis to ensure:

- Peak and motif enrichment consistent with biological function (A, C)
- High signal over background in genome-wide analyses (B)
- Reliable performance across production lots
- FOXA1 antibody (EpiCypher 13-2001) data are shown

## High quality antibodies enable low-input autoCUT&RUN experiments


**Robust profiling for multiple targets down to 10,000 nuclei**

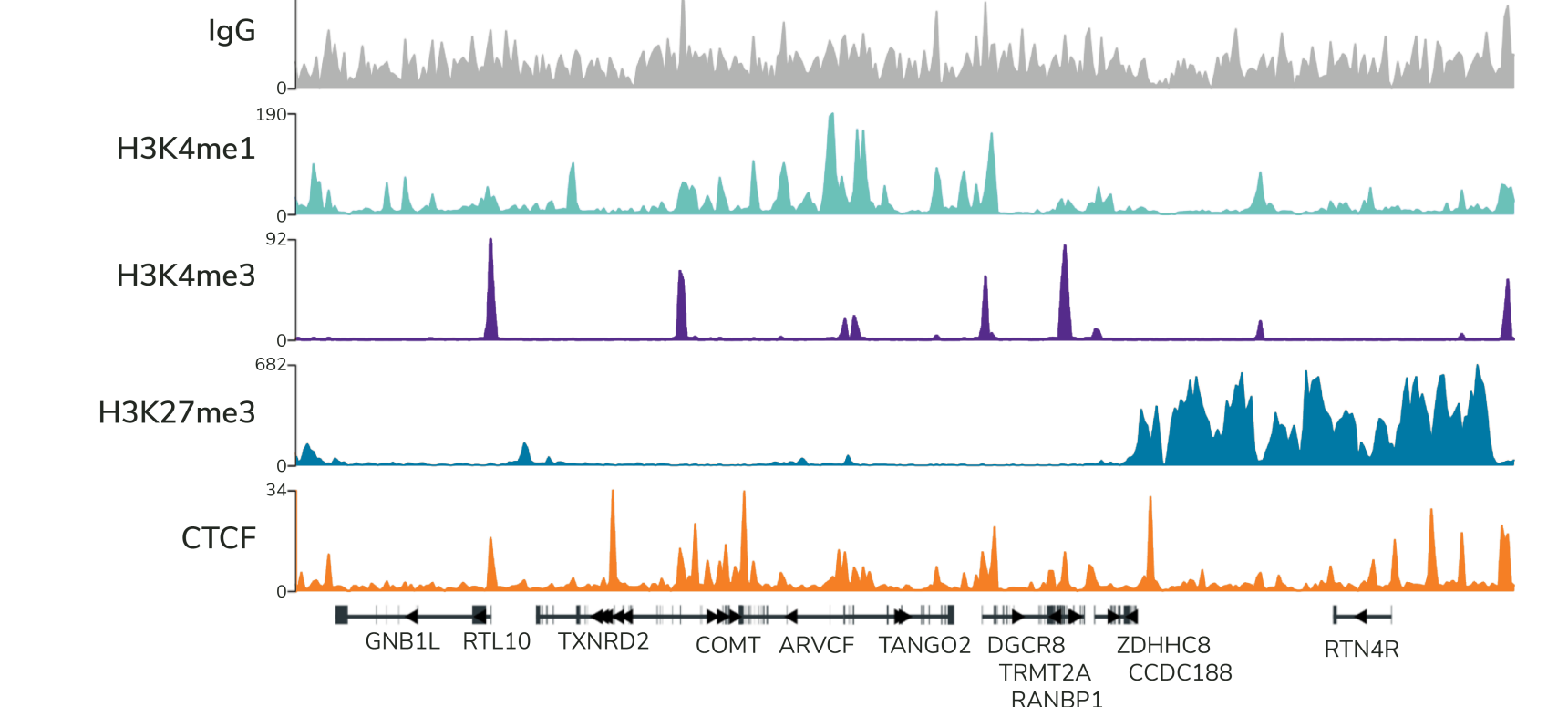

**Figure 7:** Genomic tracks were generated by autoCUT&RUN using 10,000 lightly cross-linked K562 nuclei and EpiCypher-validated antibodies to H3K4me1 (active regulatory elements), H3K4me3 (active promoters), H3K27me3 (repressed genes), and CTCF (widely studied transcription factor). See [epicypher.com/antibodies-complete-list/](http://epicypher.com/antibodies-complete-list/) for antibodies.

## Proof-of-concept: Robust profiling using 10,000 mouse immune cells

**Optimized for lightly cross-linked nuclei from mouse B cell line**

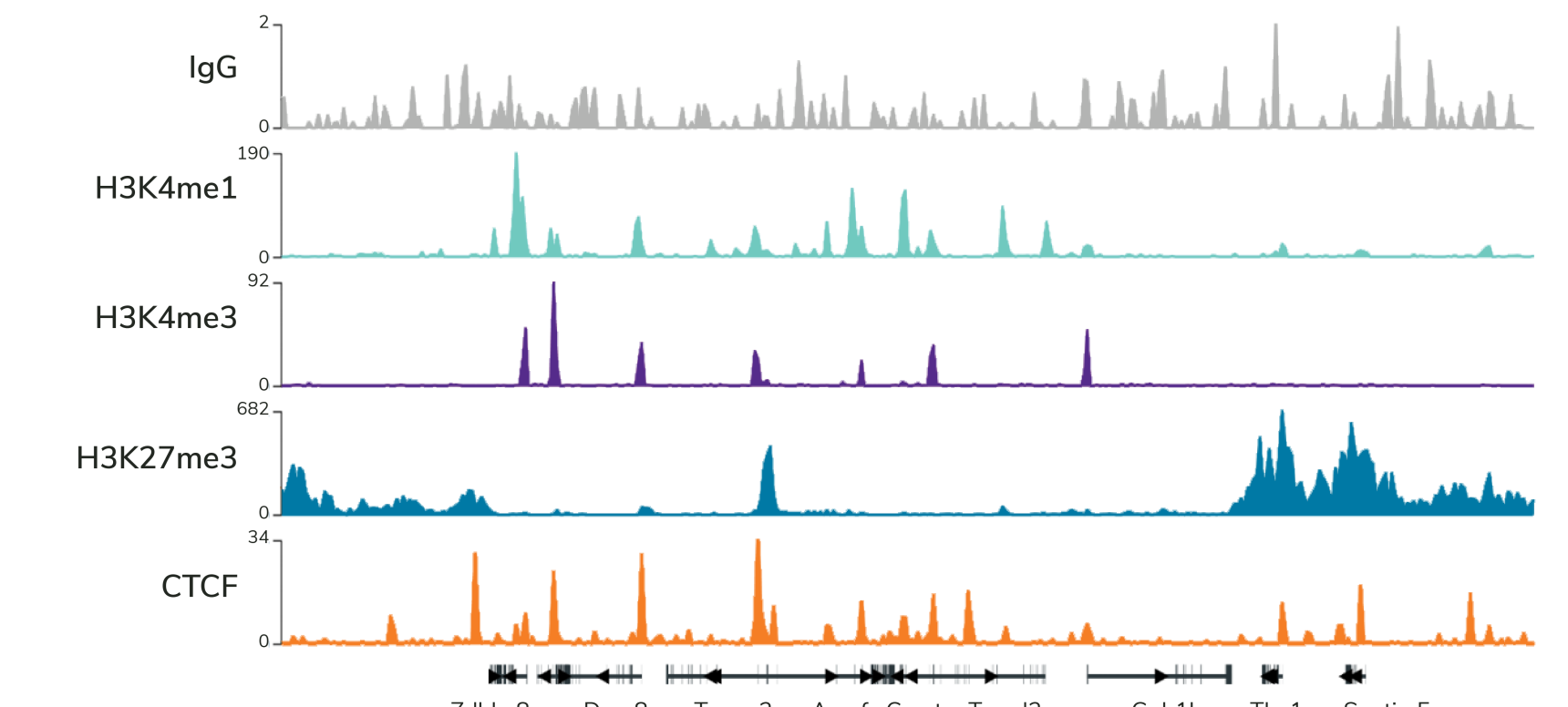

**Figure 8:** autoCUT&RUN is reliable for immune cell profiling. autoCUT&RUN was performed using 10,000 lightly cross-linked mouse B cell nuclei (provided by the ImmGen Consortium) and same antibodies as in Figure 7.


## Conclusions

- **Epigenomics is central to understanding gene regulatory processes**, but historical methods (ChIP-seq) are unreliable
- Ultra-sensitive CUTANA CUT&RUN assays are poised to dramatically change the field, improving access to high-resolution chromatin mapping
- EpiCypher developed **automated CUT&RUN for high throughput and cost-effective chromatin mapping**, ideal for studying cell & gene therapy at scale
- These efforts are bolstered by **quantitative SNAP-CUTANA Spike-in Controls and highly specific antibodies**, both of which were crucial to optimizing autoCUT&RUN for immune cell profiling

# autoCUT&RUN defines immune cell differentiation pathways for advanced cell & gene therapy research


- Collaboration with ImmGen Consortium
- Each assay only required 10,000 FACS-sorted cells
- Compatible with stimulated and cross-linked cells
- Standardized sample handling improved yields
- Cell-type specific peaks detected – useful for studying cell differentiation

## Robust data across diverse primary mouse immune cell types

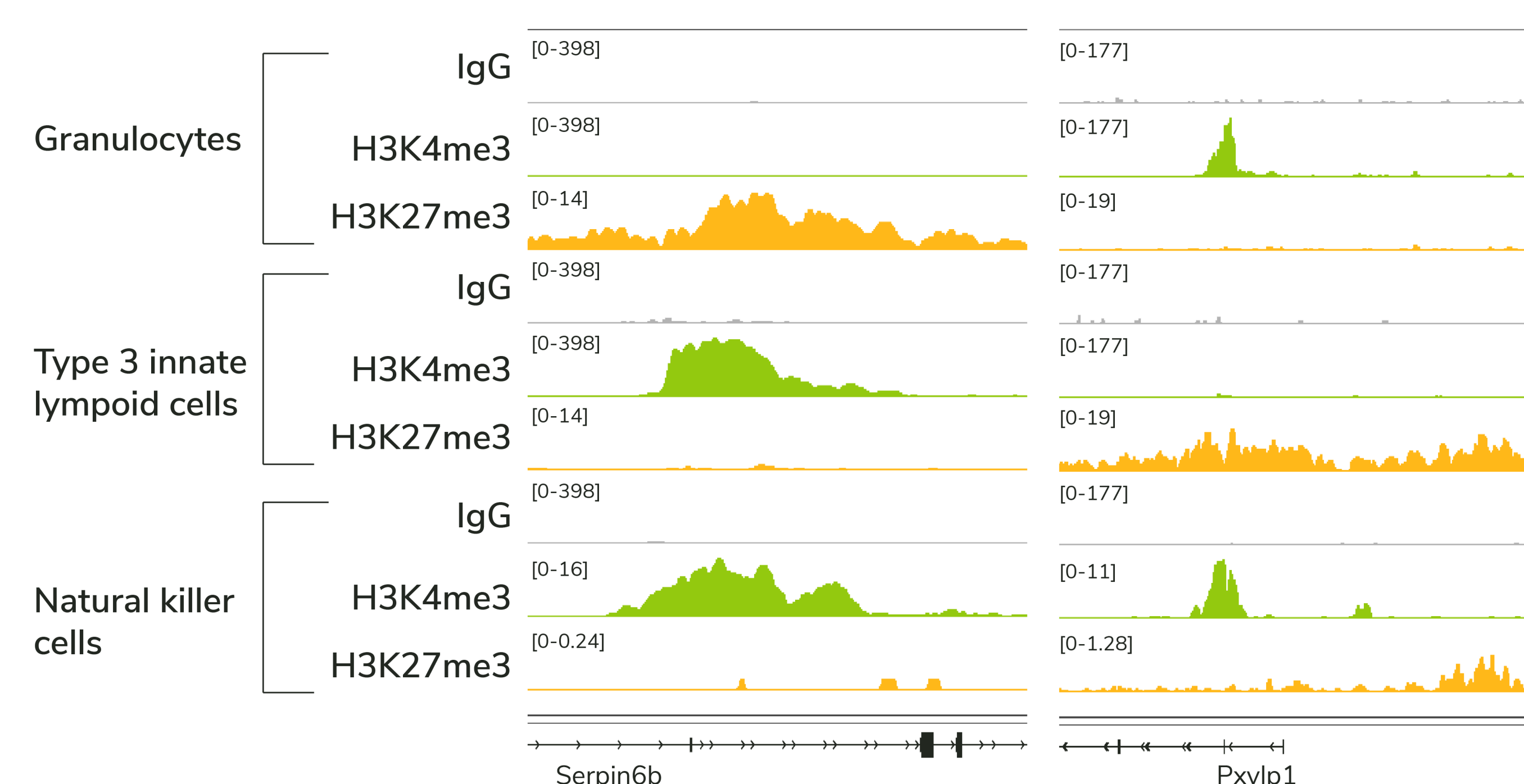

**Figure 9.** autoCUT&RUN reveals distinct H3K4me3 (marks active promoters) and H3K27me3 (denotes repressed genes) profiles across FACS-sorted primary mouse granulocytes, type 3 ILCs, and NK cells (Ly49H+), provided by ImmGen. 10,000 nuclei were used per reaction.

## In-depth studies of individual immune cell populations reveals novel biology

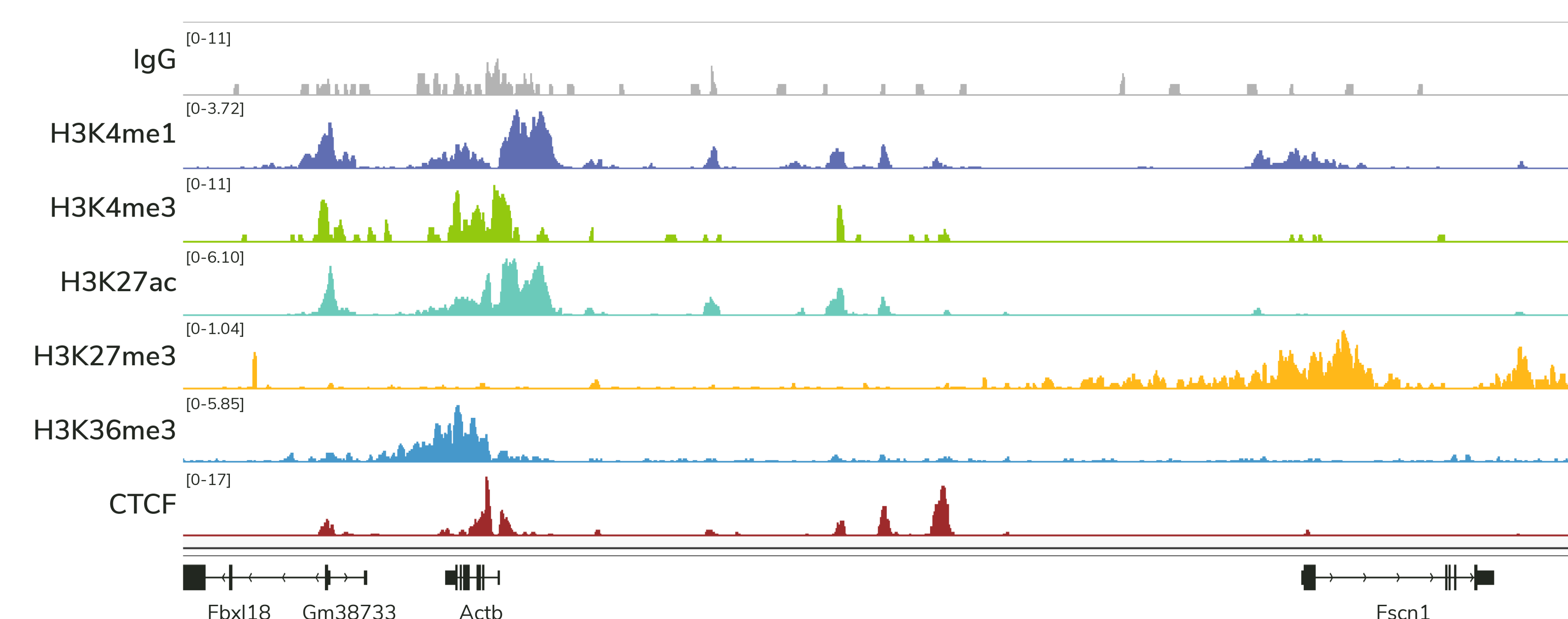

**Figure 10.** High-resolution profiling of FACS-sorted type 3 ILCs using autoCUT&RUN identifies unique genomic compartments, including active regulatory elements (H3K4me1, H3K27ac), promoters (H3K4me3), and gene bodies (H3K36me3), as well as repressed genes (H3K27me3) and transcription factor binding sites (CTCF). 10,000 nuclei (from ImmGen) were used per reaction.

## Applications of CUTANA assays

- Exhausted T cells (PMID: 35930654)
- CAR T cell expansion (PMID: 36944333)
- dCas9/Cas9 targeting (PMID: 35849129)

